
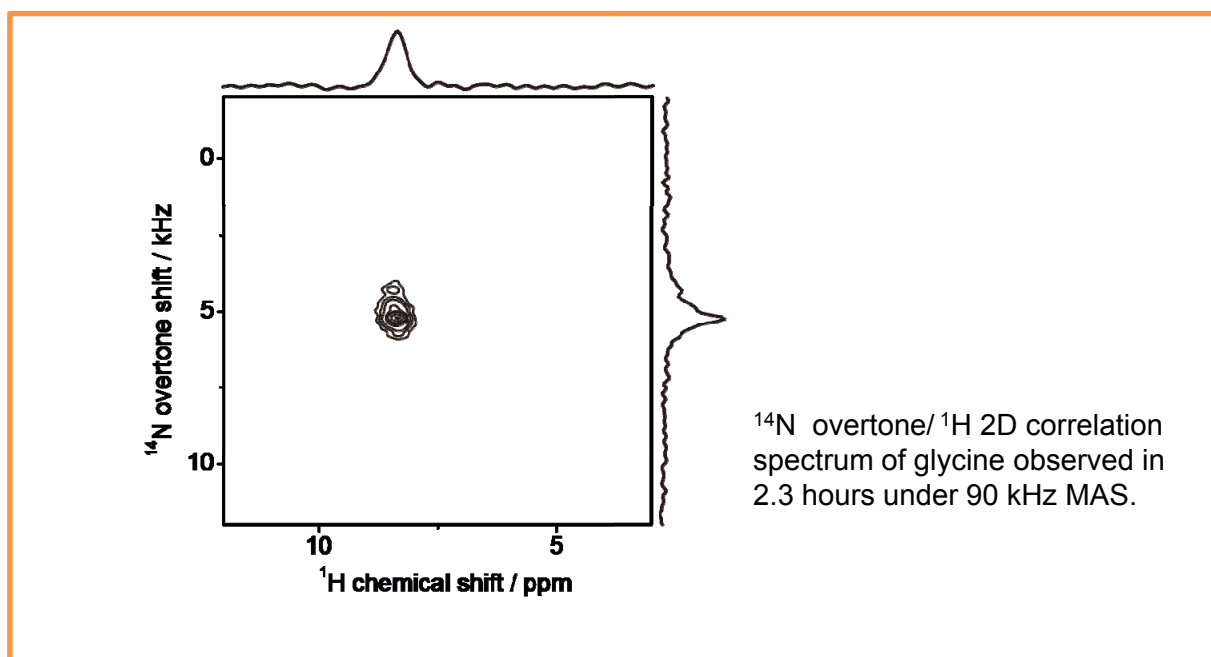


Overtone solid-state NMR spectroscopy on Nitrogen-14

Product used : Nuclear Magnetic Resonance Spectrometer (NMR)

¹⁵N NMR is widely used because of the importance of nitrogen in chemistry, materials, biology, environment etc. However, very low abundance of ¹⁵N (<0.4%) results in poor sensitivity and thus makes observation time-consuming. On the other hand, the rest of nitrogen atoms are also NMR sensitive nucleus of ¹⁴N. Despite the high abundance of ¹⁴N (>99%), it's application is rather limited due to the huge quadrupolar interactions and its spin quantum number I = 1. The introduction of very fast MAS (>70 kHz) enabled the ¹H detected ¹⁴N/¹H correlation spectroscopy (Nishiyama et al., JMR 208 (2011) 44-48). However, it demands precise magic-angle adjustment, moreover, the sensitivity is largely affected by the molecular motion. It was shown that the use of ¹⁴N overtone transition avoids these difficulties under MAS (O'Dell et al. Chem. Phys. Lett. 514 (2011) 168-173) with a cost of sensitivities. Since the overtone transition between -1 and +1 energy level is forbidden, the low sensitivity fundamentally comes from small transition moment between these levels. Here, we have developed ¹H detected ¹⁴N overtone/¹H correlation spectroscopy under ultrafast MAS conditions > 70 kHz to achieve highly sensitive ¹⁴N NMR spectroscopy of rigid solid.

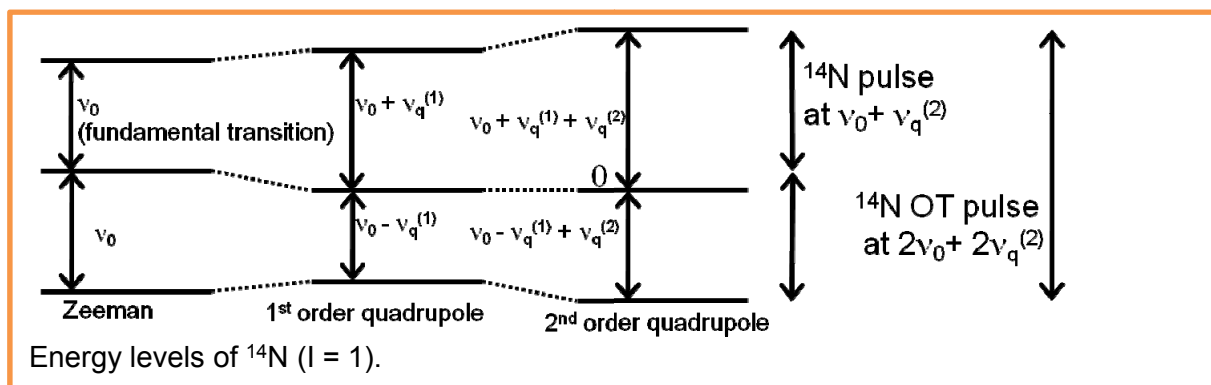

Y. Nishiyama, M. Malon, Z. Gan, Y. Endo, T. Nemoto, J. Magn. Reson. 230 (2013) 160-164.

Copyright © 2018 JEOL Ltd.

Certain products in this brochure are controlled under the "Foreign Exchange and Foreign Trade Law" of Japan in compliance with international security export control. JEOL Ltd. must provide the Japanese Government with "End-user's Statement of Assurance" and "End-use Certificate" in order to obtain the export license needed for export from Japan. If the product to be exported is in this category, the end user will be asked to fill in these certificate forms.

