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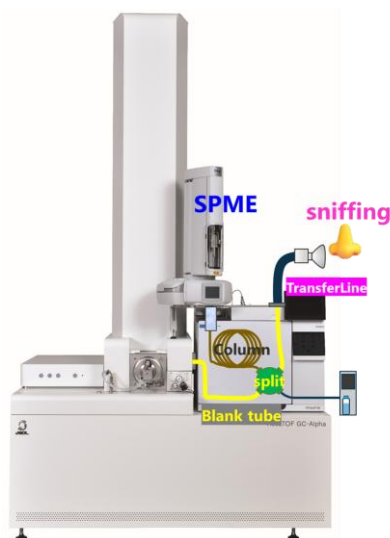
Figure 1 JMS-T2000GC with HT2850T autosampler and schematic diagram of sniffing

Table 1 Measurement condition

HS-SPME conditions		Sniffing Conditions	
Auto-sampler	HT2850T (HTA S.R.L.)	Sniffing port	OP275 (GL Sciences Inc.)
SPME fiber	DVB/CAR/PDMS 2cm (MERCK)	Transfer line temperature	200°C (constant)
Extraction	50 °C, 30min	Flow ratio (transfer line : MS)	1:1 (oven temp. 40°C), 7:3 (oven temp. 250°C)
Desorption	250 °C, 3min		
GC conditions		MS conditions	
Gas Chromatograph	8890 GC (Agilent Technologies)	Spectrometer	JMS-T2000GC (JEOL Ltd.)
Column	DB-5MS (Agilent Technologies) 30m x 0.25mm, 0.25µm	Ion source	EI/FI combination
Inlet	250 °C, Splitless	Ionization	EI(70eV, 250 °C), FI
Oven temperature	40°C(3min) → 10°C/min → 250°C(5min)	Mass range	m/z 10-800
Carrier flow	He, 131 kPa (Constant Pressure)	Analysis software	msFineAnalysis AI

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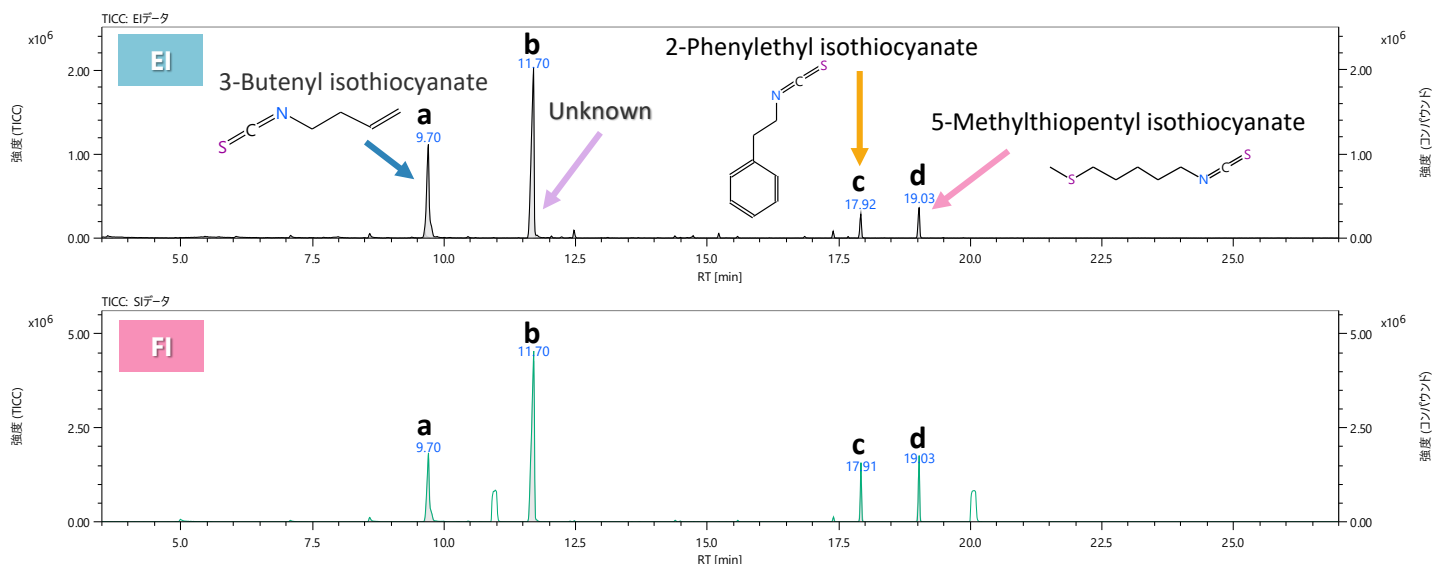
Figure 2 Measurement results of pickled Chinese cabbage: TIC chromatogram

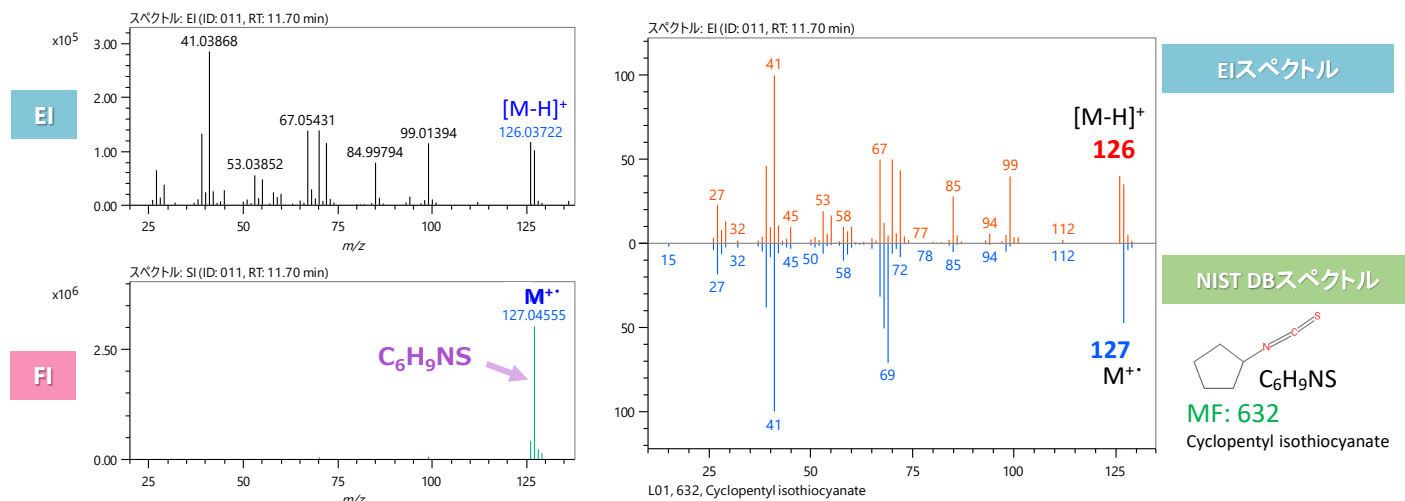
Figure 3 Comparison of the mass spectrum of compound B with that of the NIST DB search results

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Figure 4 AI structural analysis results window for compound b

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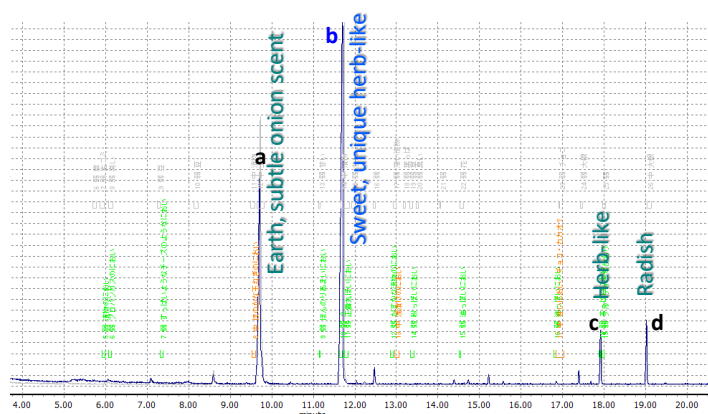
Figure 5 Integration of total ion current chromatogram (TIC) and odor information in GC-O analysis results

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