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Figure 1 JMS-T2000GC AccuTOF™ GC-Alpha 2.0

Table 1 GCxGC-MS measurement condition

GCxGC		MS	
Column	1st: BPX5, 30 m × 0.25 mm, 0.25 μm, 2nd: Rxi17-SilMS, 3.4 m × 0.15 mm, 0.15 μm	Ion Source	El standard ion source / CI ion source
Inlet	250 °C, Split 10:1	Ionization	El+: 70 eV, 300 μA, CI+: 200 eV, 300 μA, Reagent Gas: isobutane
Oven	80 °C (2 min) → 5 °C/min → 325 °C (9 min)	Acq. Speed	El: 50 Hz, CI: 25 Hz
Carrier flow	He, 1.5 mL/min (Constant Flow)	m/z Range	El: m/z 33 - 800, CI: m/z 100 - 800
Modulation period	6 sec		

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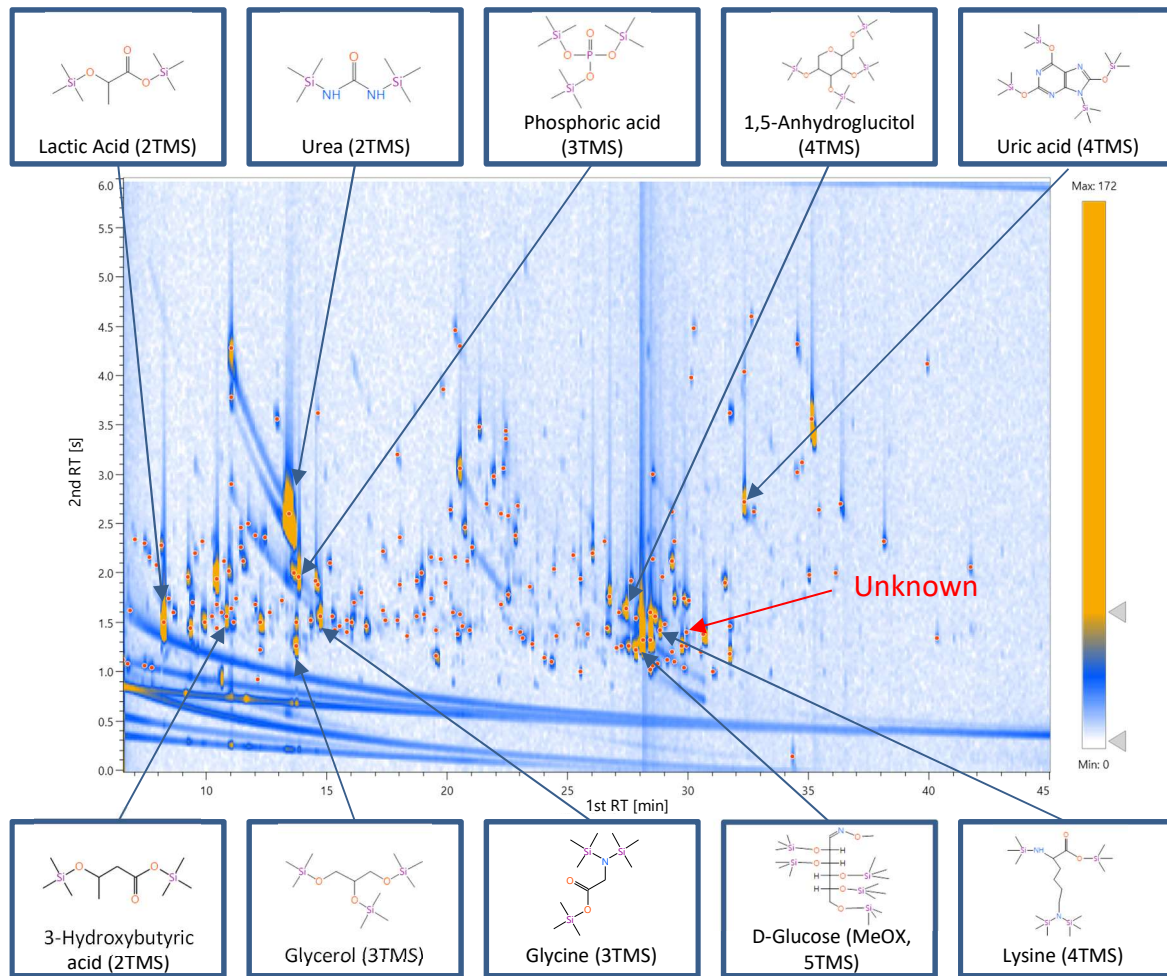
Figure 1 GCxGC TIC of metabolites in NIST SRM 1950

Table 2 Water-soluble metabolites listed in the NIST SRM 1950 COA and detected in this measurement

#	Derivatized Compound Name	Original Compound	Certified Category	#	Derivatized Compound Name	Original Compound	Certified Category
1	Alanine_2TMS	Alanine	Certified Values	14	Tyrosine_2TMS	Tyrosine	Certified Values
2	Glycine_3TMS	Glycine	Certified Values	15	Tyrosine_3TMS	Tyrosine	Certified Values
3	Isoleucine_1TMS	Isoleucine	Certified Values	16	Valine_TMS	Valine	Certified Values
4	Isoleucine_2TMS	Isoleucine	Certified Values	17	Valine_2TMS	Valine	Certified Values
5	Leucine_1TMS	Leucine	Certified Values	18	Glucose_MOX_5TMS_PK1	Glucose	Certified Values
6	Leucine_2TMS	Leucine	Certified Values	19	Glucose_MOX_5TMS_PK2	Glucose	Certified Values
7	Lysine_3TMS	Lysine	Certified Values	20	Urea_2TMS	Urea	Certified Values
8	Lysine_4TMS	Lysine	Certified Values	21	Uric acid_4TMS	Uric Acid	Certified Values
9	Methionine_2TMS	Methionine	Certified Values	22	Phenylalanine_TMS	Phenylalanine	Non-Certified Values
10	Proline_1TMS	Proline	Certified Values	23	Phenylalanine_2TMS	Phenylalanine	Non-Certified Values
11	Proline_2TMS	Proline	Certified Values	24	Threonine_2TMS	Threonine	Non-Certified Values
12	Serine_2TMS	Serine	Certified Values	25	Threonine_3TMS	Threonine	Non-Certified Values
13	Serine_3TMS	Serine	Certified Values				

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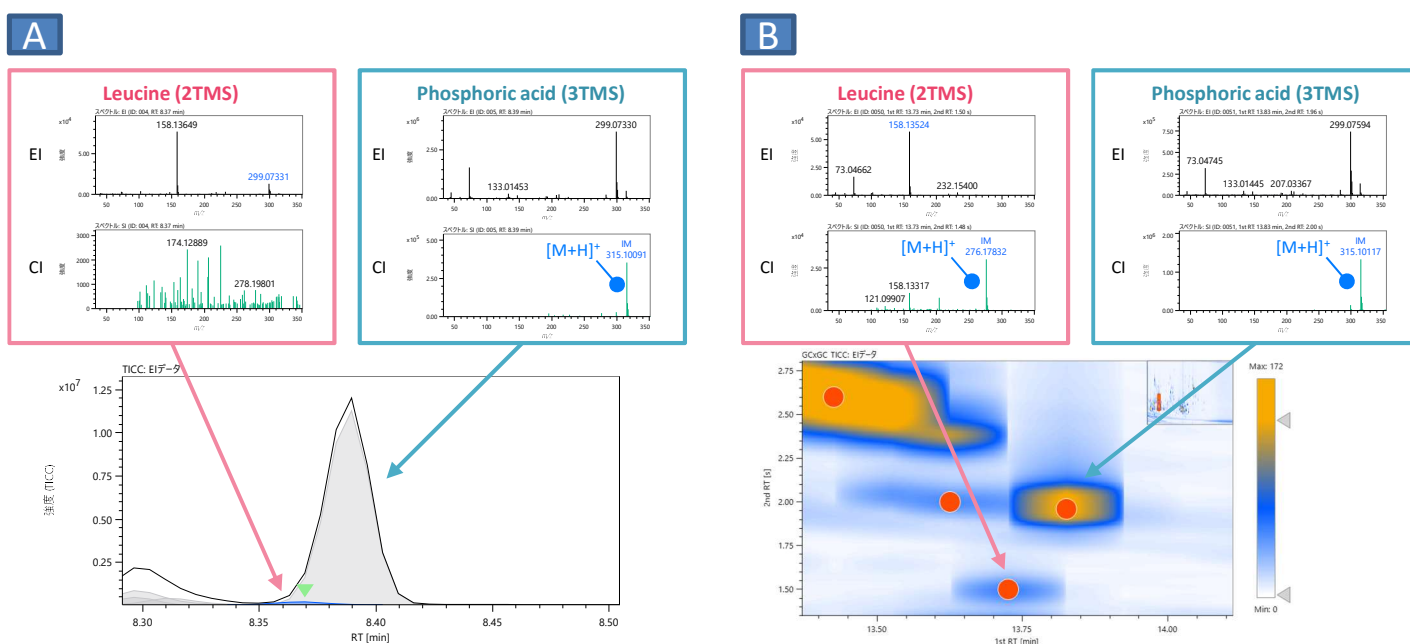
Figure 2 TICC and mass spectra of Leucine (2TMS) and Phosphoric acid (3TMS)  
A: One dimensional GC-MS result, B: GCxGC-MS result.

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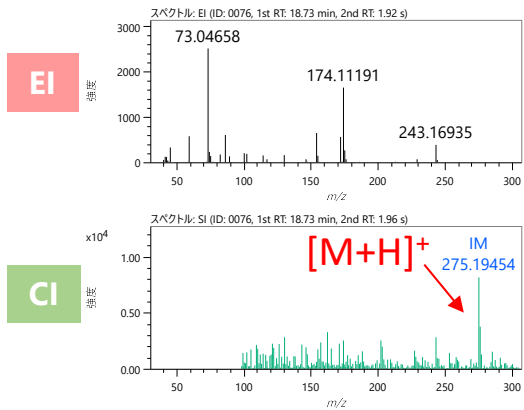
Figure 3 Mass spectra of unknown compound

Table 3 Integrated analysis result of unknown compound

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Structural formula of Acetamide, 2-amino-N-cyclopropyl-, 2TMS:

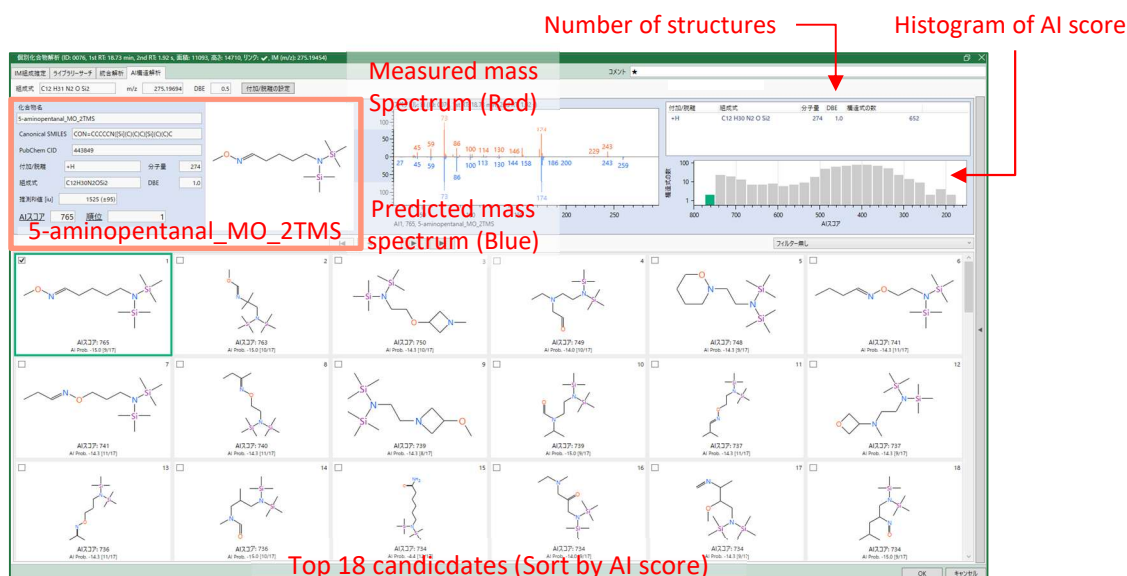
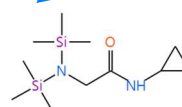
Figure 4 AI structure analysis result of unknown compound

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